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
PURELAB[®] Chorus I

Operator Manual


MANU39998 VER 02 3/17

Life Science / Analytical Research / General Science


WELCOME

Thank you for purchasing the PURELAB® Chorus 1. Please read the following Health and Safety notice before proceeding.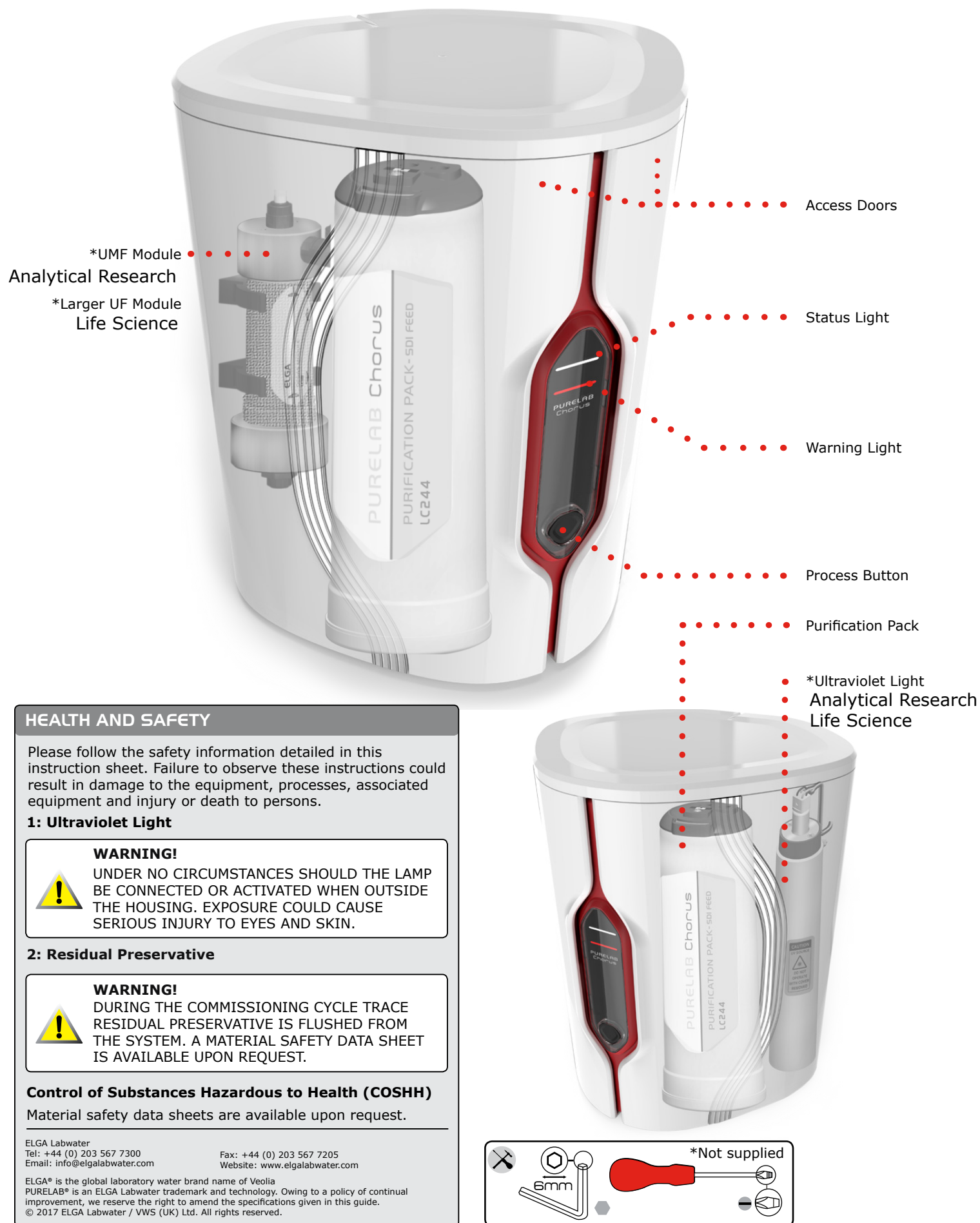


SYSTEM CONFIGURATION Life Science / Analytical Research / General Science


Example one: 1 x PURELAB® Chorus 3 + Reservoir + Chorus 1 + 1 X Dispense.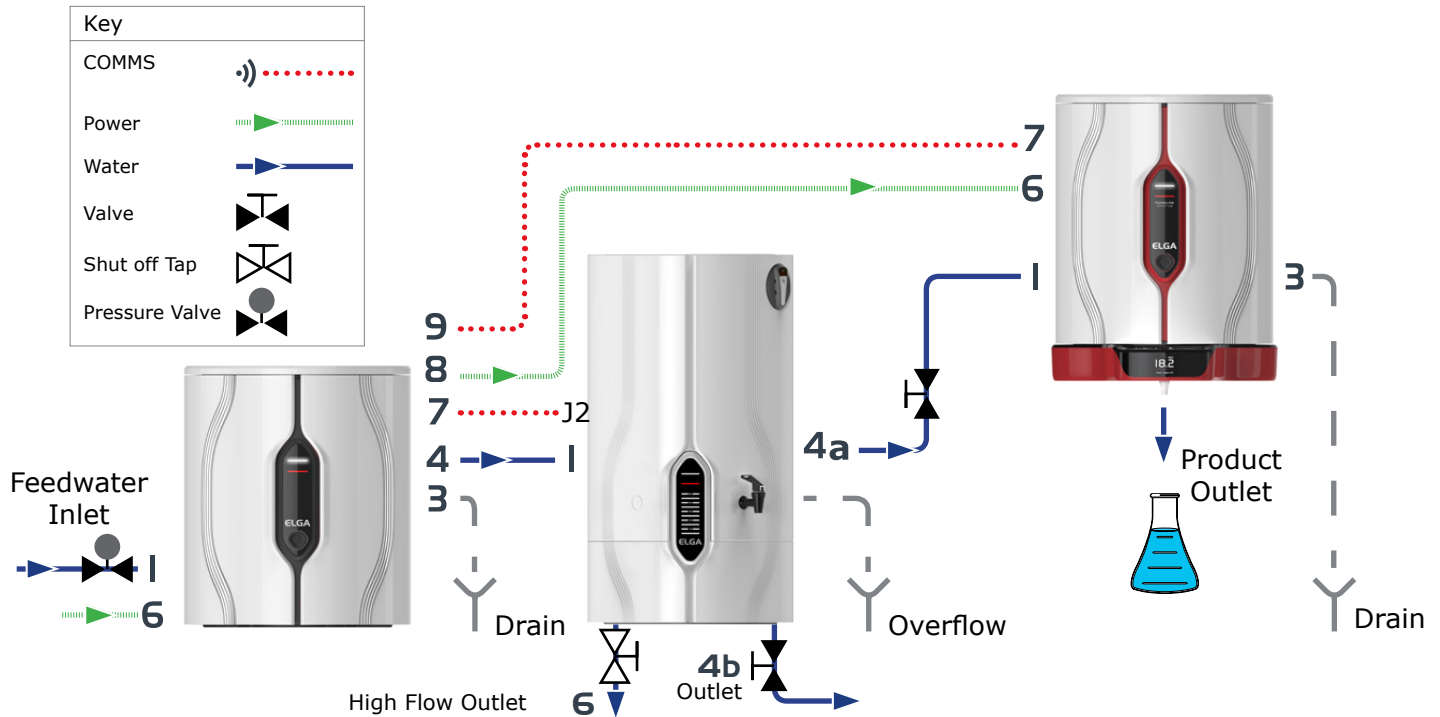

Example two: 1 x PURELAB® Chorus 1 + 2 X Dispenser.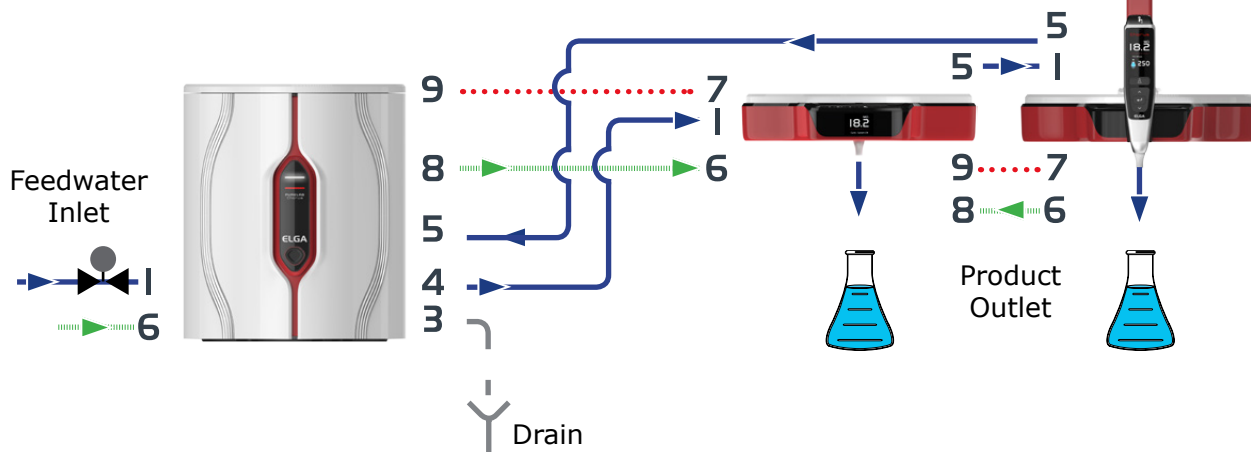


Example three: 1 x PURELAB® Chorus 1 + 3 X Dispenser.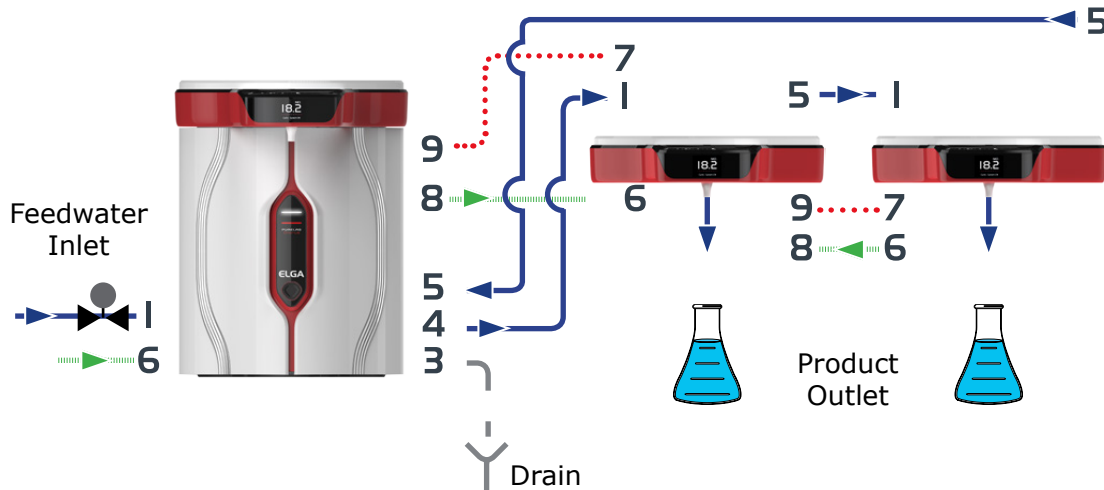

Further advice on configuration connections.

Ia MODULAR STACKING BENCH MOUNTED RESERVOIR ON TOP


PC1LSCXM2 / PC1ANRXM2 / PC1GSCM2

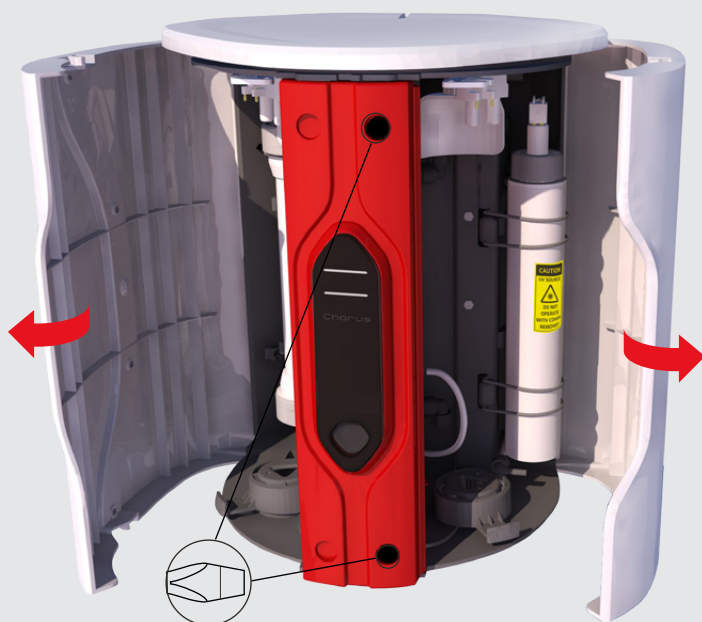a. Open doors and undo quarter turn fasteners.

b. Remove front panel and undo lid bolt with an allen key. (6mm AF X FXBOYR301716)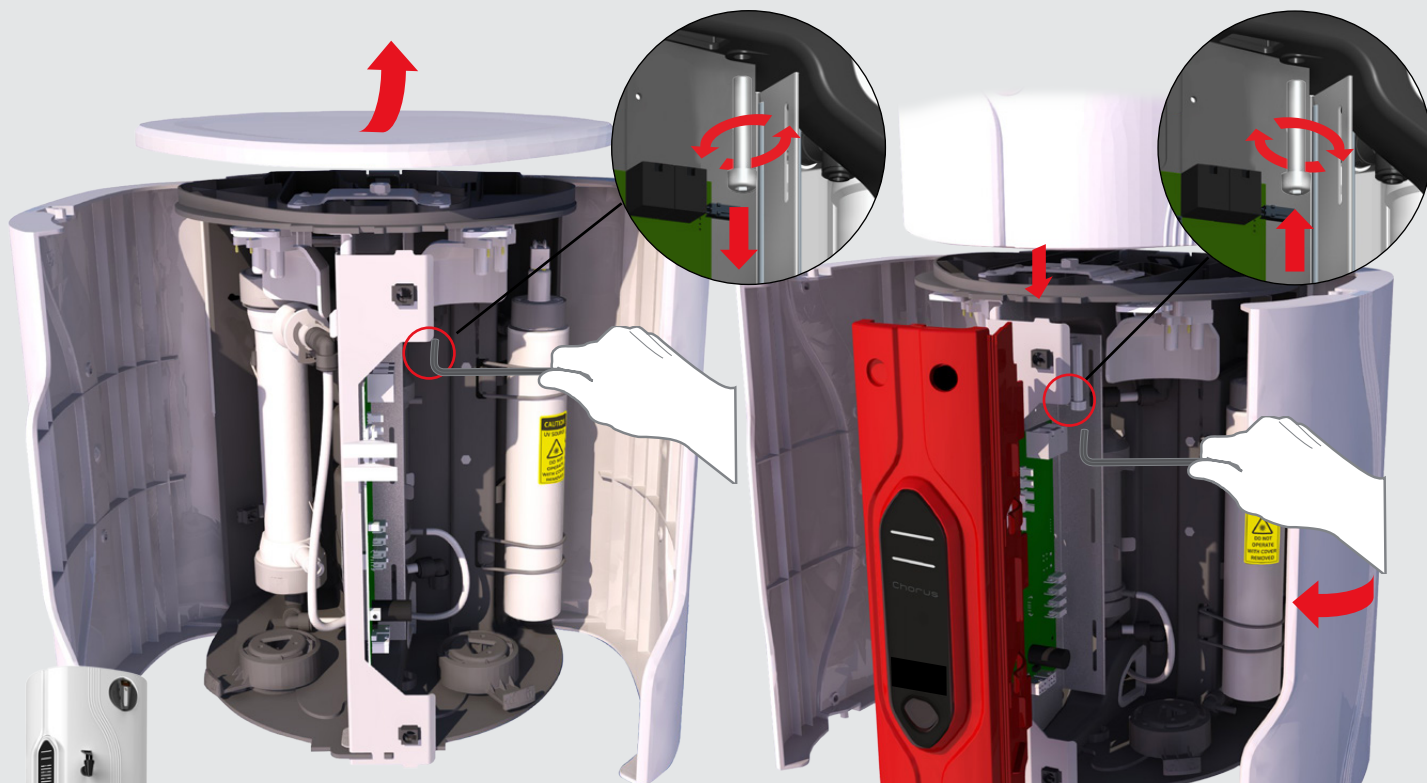



c. Remove bolt and lift the lid.

d. Place the reservoir on top, lining up with the three points. Screw in allen head bolt, before replacing front panel.


NB. Please refer to the dispenser Quick Start Guide to combine a dispenser.

**WARNING!**

ONLY PLACE ONE CHORUS SYSTEM ON TOP OR BELOW THE RESERVOIR.

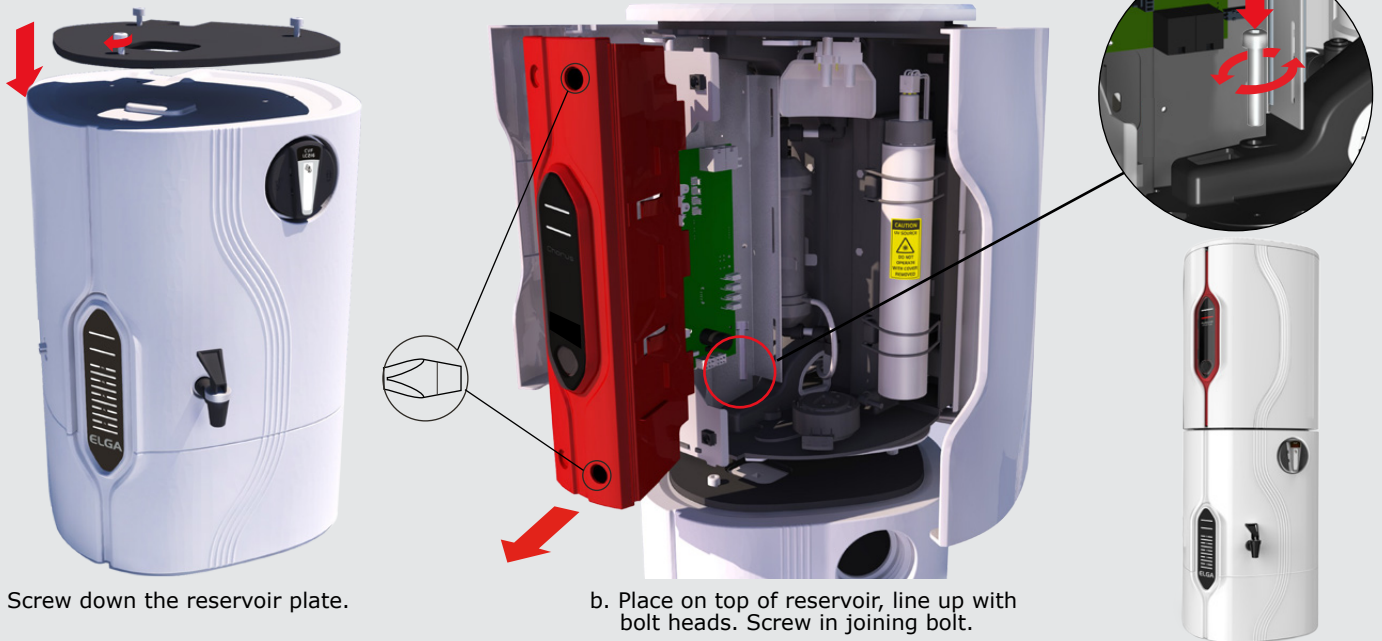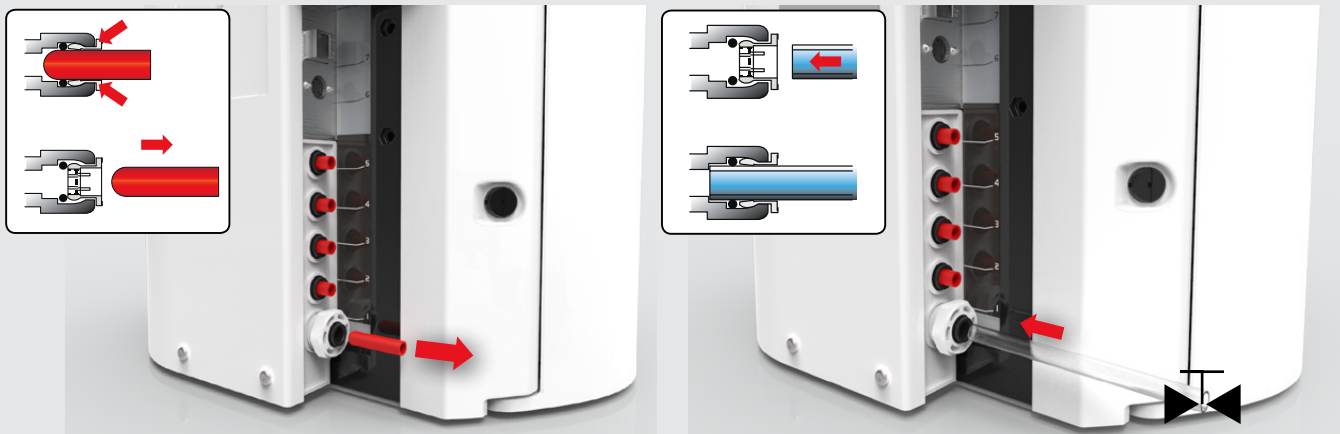
1b MODULAR STACKING RESERVOIR UNDERNEATH**2 CONNECT THE FEED WATER (PORT NO.1)****3 CONNECT THE DRAIN (PORT NO.3)****CAUTION!**

Ensure drain pipe is not in waste water.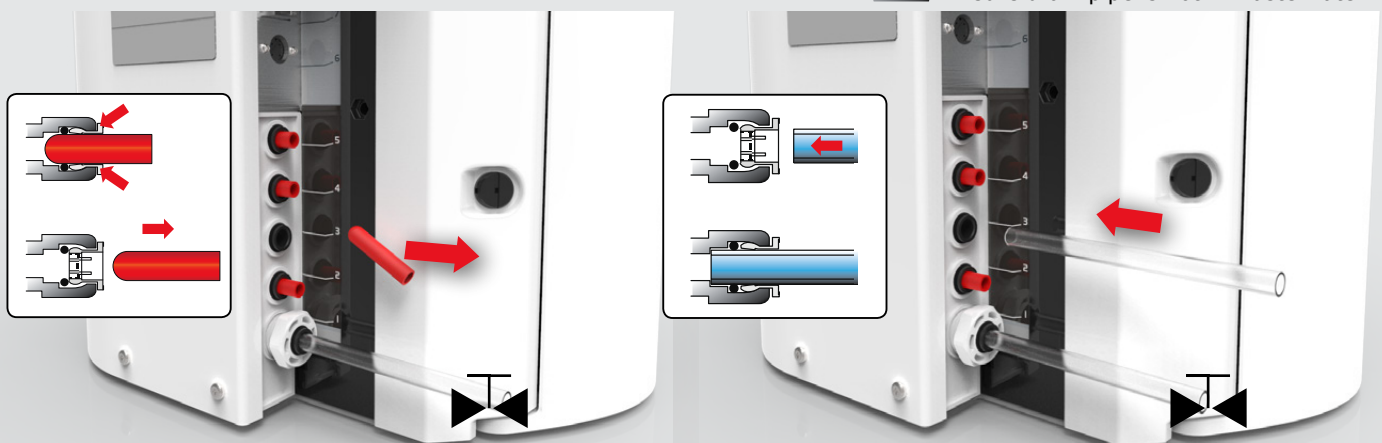


4 RETURN DISPENSER

NB: Please refer to dispenser manual for more information.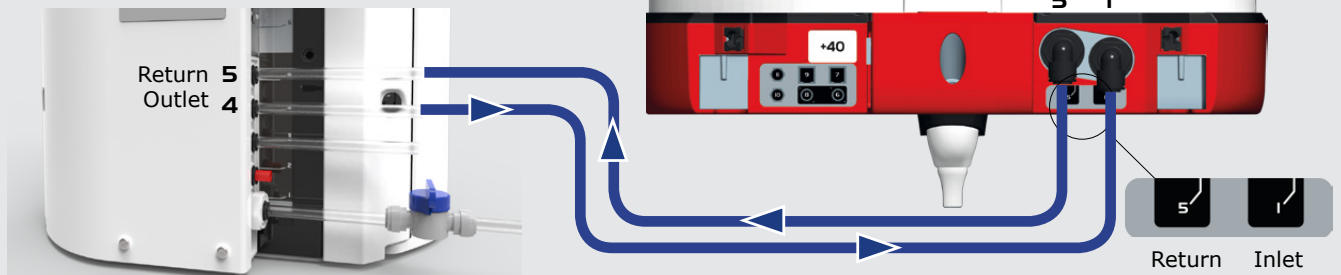

5 CONNECT THE PURIFIED WATER OUTLET (PORT NO.4)

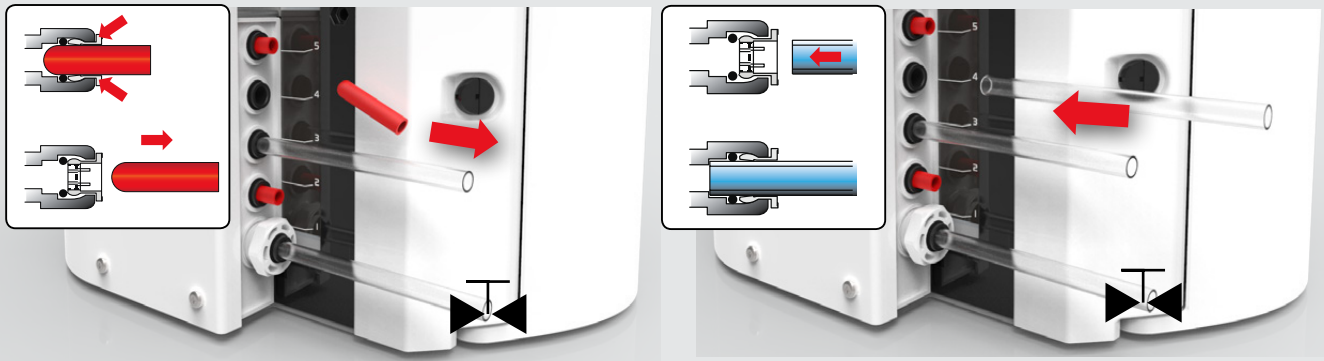a. Remove the transit plugs from the water outlet connection.

b. Using the tube supplied, firmly push one end securely into the water connector.

6 COMMS CONNECTIONS

WARNING!

ONLY USE COMMS CABLE SUPPLIED BY ELGA.

NB. Please refer to the dispenser Quick Start Guide to combine a dispenser.

7 FOLLOW THE INSTRUCTIONS

a. Open right hand door.


b. Insert Purification Pack and turn handle.

c. Make sure it's secure.

8 CONNECT THE POWER (NO.6 PORT)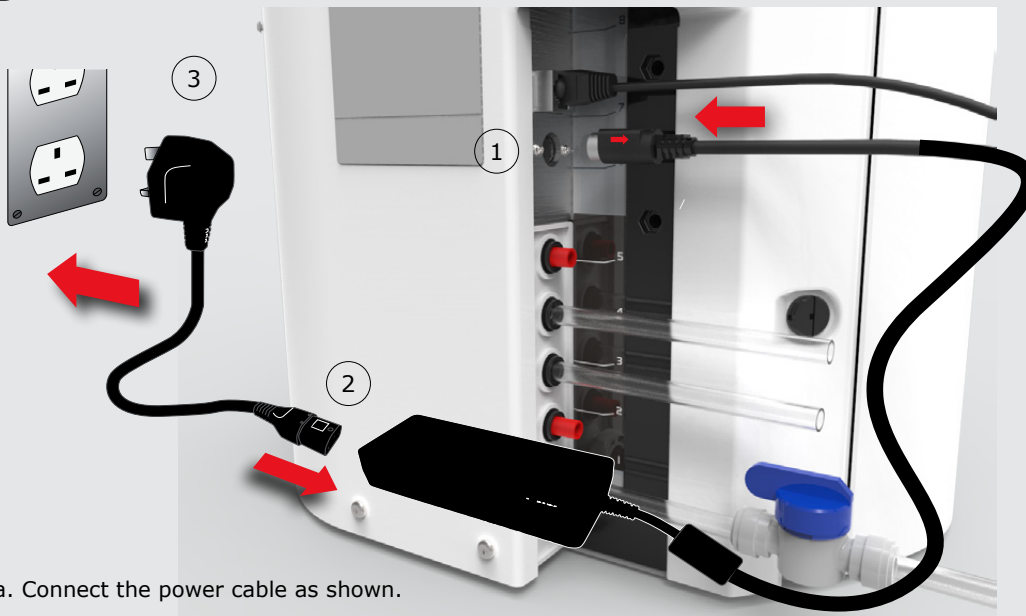

a. Connect the power cable as shown.

b. Only use the power supply supplied with the Chorus.
The power supply must be earthed.

c. **PRESS ONCE**
to start Commission mode.
Bottom red light will be
flashing until cycle has ended.

Approx 80 minutes

WARNING!

ONLY USE THE APPLIANCE COUPLER (MAINS LEAD) AND POWER SUPPLY PROVIDED. IF THE EQUIPMENT IS USED IN A MANNER NOT SPECIFIED BY ELGA, THE PROTECTION PROVIDED BY THE EQUIPMENT MAY BE IMPAIRED. POSITION THE POWER SUPPLY SO THAT IT CANNOT COME INTO CONTACT WITH WATER.

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PURELAB®
Chorus

Table of contents

1.	Introduction	Page 1
1.1	Health & Safety	Page 1
1.2	Product Range	Page 1
1.3	Use of this Manual	Page 1
1.4	Installation	Page 1
1.5	System Configuration	Page 1
1.6	Environment	Page 1
1.7	Commissioning	Page 1
1.8	Customer Support	Page 1
2.	Your Guide to Chorus	Page 2
2.1	PURELAB® Chorus Part No. PC1LSCXM2 - Life Science	Page 2
2.2	PURELAB® Chorus Part No. PC1ANRXM2 - Analytical Research	Page 3
2.3	PURELAB® Chorus Part No. PC1GSCXM2 - General Science	Page 4
3.	Maintenance	Page 5
3.1	Cleaning the inlet filter assembly	Page 5
3.2	Replacing the Purification pack	Page 6 - 7
3.3	Replacing the Ultraviolet lamp	Page 8 - 9
3.4	Replacing the Ultra filtration module	Page 10 - 11
3.5	Replacing the Ultra mirco filtration module	Page 12 - 13
3.6	Sanitisation Procedure	Page 14
4.	Self Help Guidelines	Page 14
5.	Health & Safety notes	Page 15
5.1	Environment	Page 15
5.2	Electricity	Page 15
5.3	Residual Biostat	Page 15
5.4	Ultraviolet lamp	Page 15
5.5	Control Of Substances Hazardous to Health (COSHH)	Page 15
6.	Technical Specifications	Page 16 - 17
7.	Customized Operation	Page 17
7.1	USB Logging	Page 17
8.	Warranty / Conditions	Page 18

Copyright Note

The information contained in this document is the property of VWS (UK) Ltd, trading as ELGA® LabWater, and is supplied without liability for errors or omissions.

No part of this document may be reproduced or used except as authorized by contract or other written permission from VWS (UK) Ltd. The copyright and all restrictions on reproducing and use apply to all media in which this information may be placed.

VWS (UK) Ltd. Pursue a policy of continual product improvement and reserve the right to alter without notice the specification, design, price or conditions of supply of any product or service.

© VWS (UK) Ltd. 2017 - All rights reserved.
ELGA® and PURELAB® Are registered trademarks of VWS (UK) Ltd.

Copyright Note

Effersan™ is a registered trademark of Activon Inc.

INTRODUCTION

1.1 Health & Safety

Please ensure you read the Health & Safety notes in Section 5.

1.2 Product Range

This Operator manual has been prepared for PURELAB® Chorus product models:

- PURELAB® Chorus Part No. PC1LSCXM2 - Life Science
- PURELAB® Chorus Part No. PC1ANRXM2 - Analytical Research
- PURELAB® Chorus Part No. PC1GSCXM2 - General Science

1.3 Use of this Manual

This manual guides you through the operation and maintenance of the Chorus allowing you to obtain a guaranteed supply of purified water to meet your requirements.

1.4 Installation

Chorus is supplied with a Quick Start Guide and Quick Reference Guide.

Quick Start Guide shows you how to install and commission the Chorus.
Quick Reference Guide shows you how perform day to day functions.

1.5 System Configuration

Maximum system size 1 X Chorus 1 + 4 X Chorus 2 or 3 + Reservoir + 4 X Dispenser.
See Quick Start Guide for typical configurations.

1.6 Environment

The Chorus unit should be installed on a flat, level surface, in a clean, dry environment.
The unit can also be wall mounted using the purpose designed wall mounting kit against a vertical wall capable of supporting the combined weight of the unit and mounting kit.

WARNING!

NOT TO BE USED IN FUME EXTRACTION ENCLOSURES.

1.7 Commissioning

Chorus is supplied in a preset commissioning mode which must be completed before purified water can be dispensed. Refer to Self Help Guidelines, section 4 Process Button.


1.8 Customer Support

If you need help with your Chorus, please call your local ELGA representative. For the address of the nearest ELGA LabWater Sales and Service office visit the country list on our website.
www.elgalabwater.com or contact ELGA Lab-Water at: Email: techsupport@elgalabwater.com


2 YOUR GUIDE TO THE CHORUS
2.1 Life Science - PC1LSCXM2

Use a slotted screw driver to release quarter turn slots on Chorus.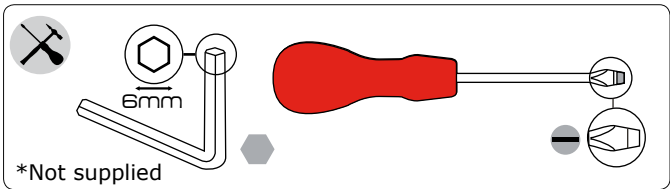


- 9. COMMS 2
- 8. Power In 2
- 7. COMMS 1
- 6. Power In 1
- 5. Return
- 4. Product Outlet
- 3. Drain
- 1. Water Inlet

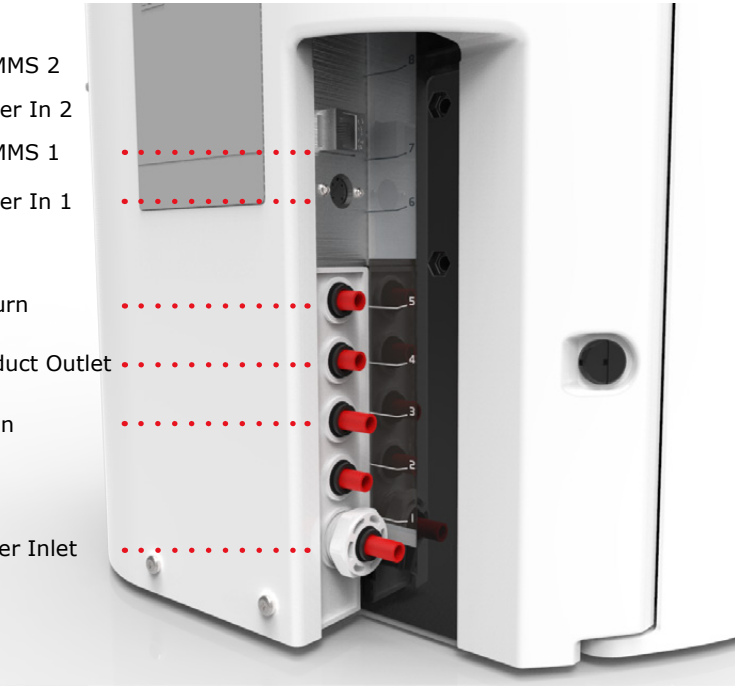2.2 Analytical Research - PC1ANRXM2

Use a slotted screw driver to release quarter turn slots on Chorus.

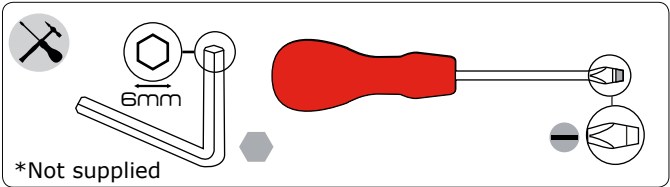
1. Water Inlet
2. Drain
3. Product Outlet
4. Return
5. COMMS 1
6. Power In 1
7. COMMS 2
8. Power In 2
9. COMMS 2


2.3 General Science - PC1GSCXM2

 Use a slotted screw driver to release quarter turn slots on Chorus.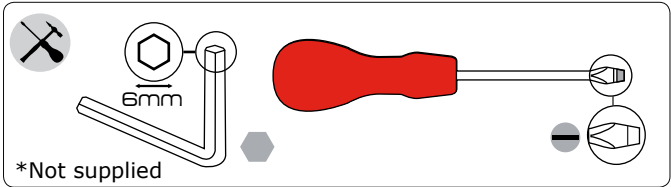

- 9. COMMS 2
- 8. Power In 2
- 7. COMMS 1
- 6. Power In 1
- 5. Return
- 4. Product Outlet
- 3. Drain
- 1. Water Inlet

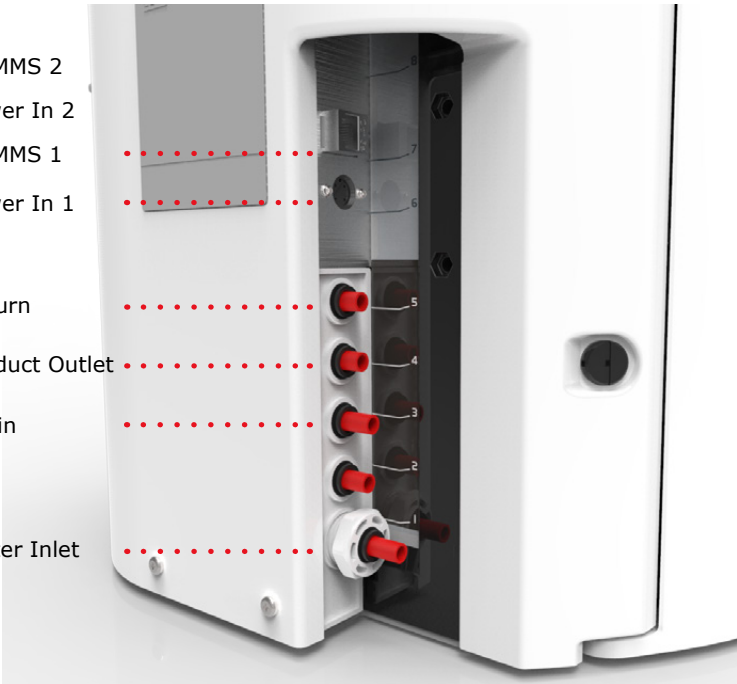3 MAINTENANCE

WARNING!


An approved supplier or distributor should carry out any maintenance work not included in this manual.

Note


Note: Disposal of all end-of-life consumable items should be in accordance with local statutory regulations.

3.1 Cleaning the Inlet Filter Assembly

Follow the instructions below.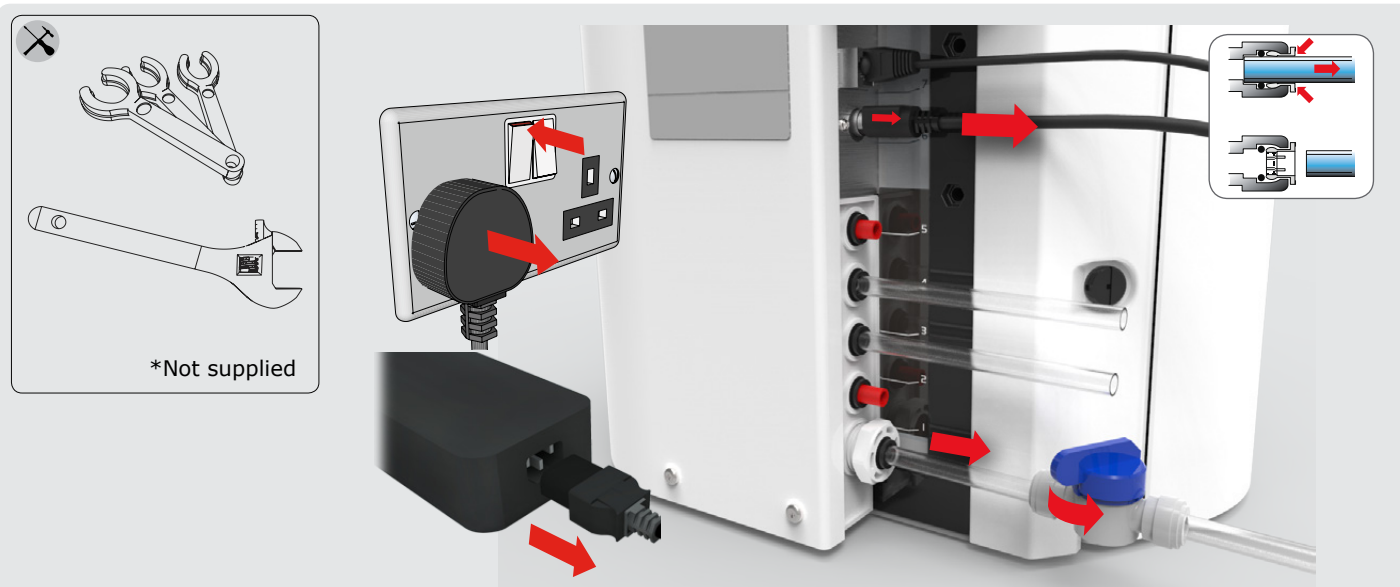

1. Isolate the electrical supply.
2. Turn off water supply, remove inlet pipe.

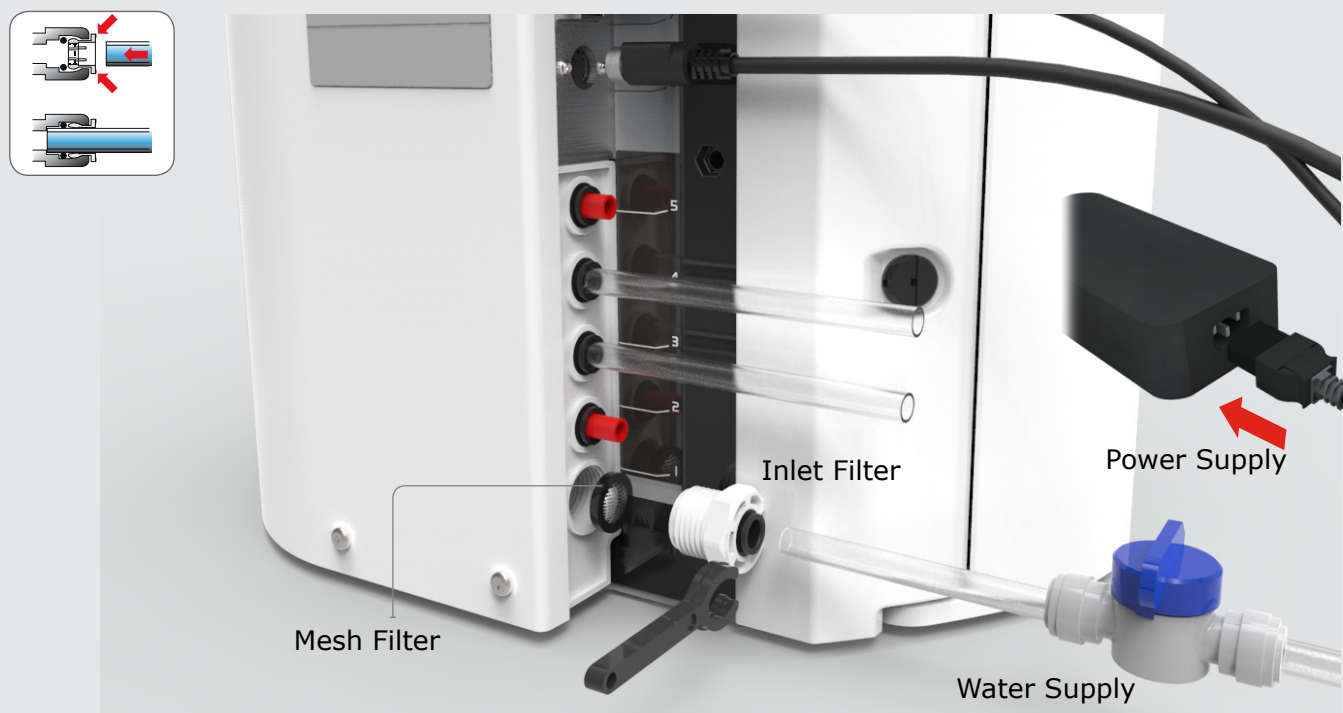3. Unscrew the inlet filter and remove the mesh filter. Replace or clean and re-assemble.
4. Reconnect the water and power supply.

3.2 Replacing the Purification Pack

Follow the instructions below.

PC1LSCXM2 / PC1ANRXM2 / PC1GSCM2

1. "Pure Sure Alarm active"
Press and hold accept to enter main menu.

2. Scroll and press accept
"Change Purification Pack(s)".

3. "Select Module" Press accept.


4. Pull down and tilt towards to remove.

5. Remove and dispose of pack 1.

6. Fit Purification Pack 2 into position 1.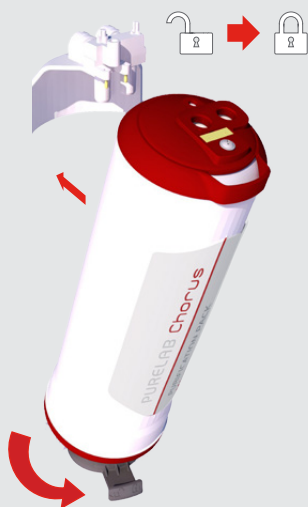

7. Insert the base first and rotate the handle.

8. Unpack New Purification Pack and remove the red transit plugs before installing.

9. Insert into position 2.

10. Insert the base first and rotate the handle. Press accept to continue.

11. Place a 5L container under the nozzle and Press dispense button to continue.

12. The system will rinse for up to 1 hour. Follow on screen instructions.

3.3 Replacing the Ultraviolet Lamp LC210

Follow the instructions below.

PC1LSCXM2 / PC1ANRXM2

Note.

The Ultraviolet (UV) lamp should be replaced when:



- UV lamp failure occurs
- UV lamp efficiency decrease affecting water purity.

**WARNING**



UV-C Radiation is harmful to the eyes and skin. UV Lamp should be operated in the chamber. It is strongly recommended that during the handling of the UV lamp cut-resistant gloves are worn.

Mercury Hazard, do not break. The lamp contains small amount of mercury. May cause redness or irritation as a result of contact with skin and / or eyes.

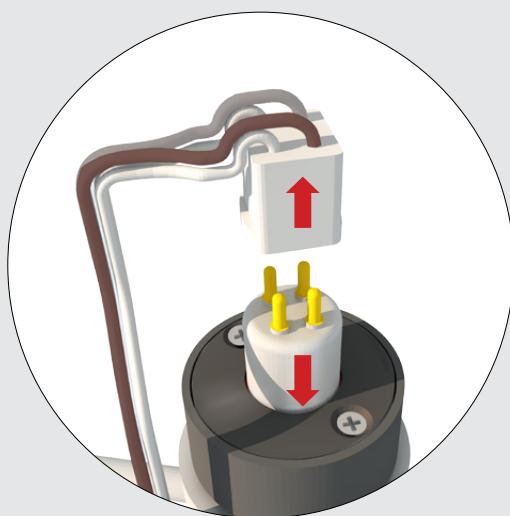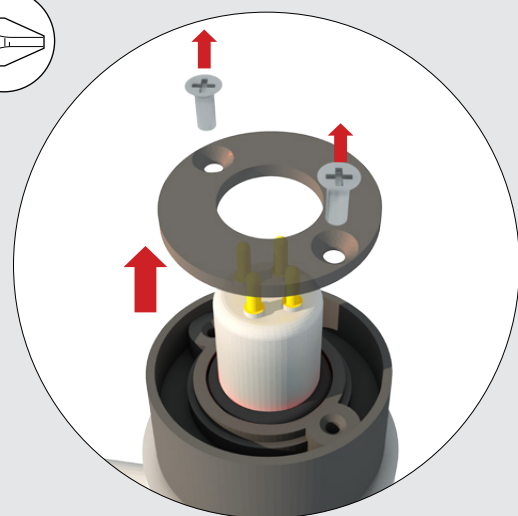**1.** Process button off, open right door.**2.** Isolate the power supply and turn off the water.**3.** Unclip UV Lamp LC210 from Chorus.**4.** Disconnect the lamp connector fitted to the top of the UV lamp.**5.** Loosen the 2 securing screws at the top of the UV housing.**CAUTION!**

Hold on to the UV lamp carefully to ensure it doesn't break.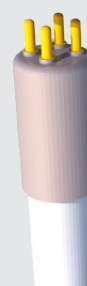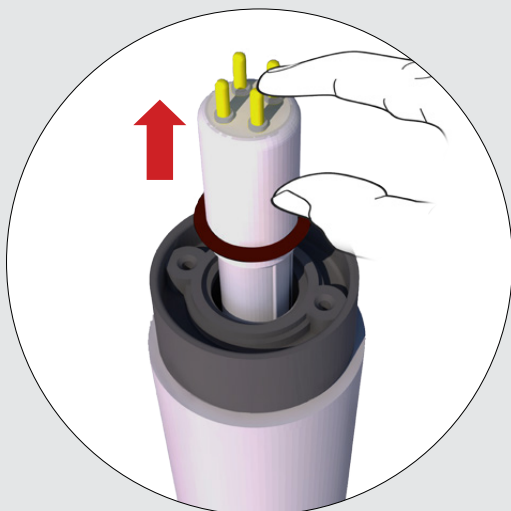

- 6.** Lift and remove the UV lamp from the housing. Discard the used UV lamp accordance with local regulations.

(Re-use the red O ring on the new UV lamp)


CAUTION!



Take care not to touch the surface of the glass. Ideally wear gloves, handle with soft cloth and wipe the surface with alcohol before fitting into the UV lamp housing.

- 7.** Unpack the new UV Lamp.

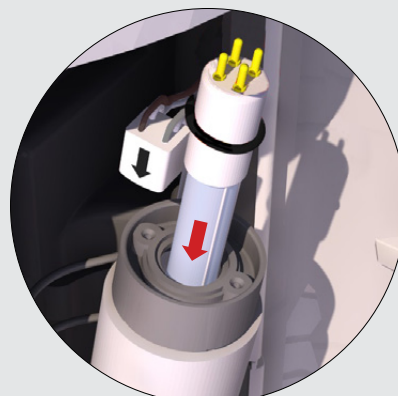
- 8.** Slide the new UV Lamp into the housing.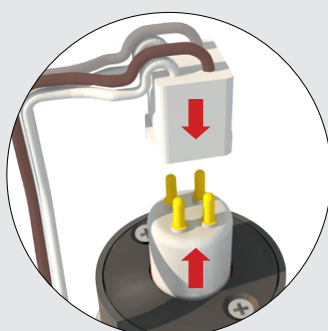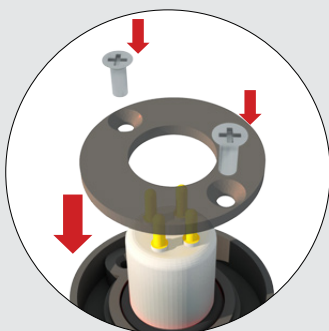

- 9.** Insert the cap and tighten the screws.

- 10.** Connect the lamp connector to the end of the UV.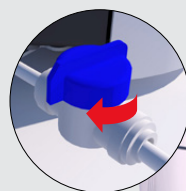

- 11.** Turn on the water supply.




- 12.** Turn on the power supply.

3.4 Replacing the Ultra Filtration Module LC151

Follow the instructions below.

PC1LSCXM2

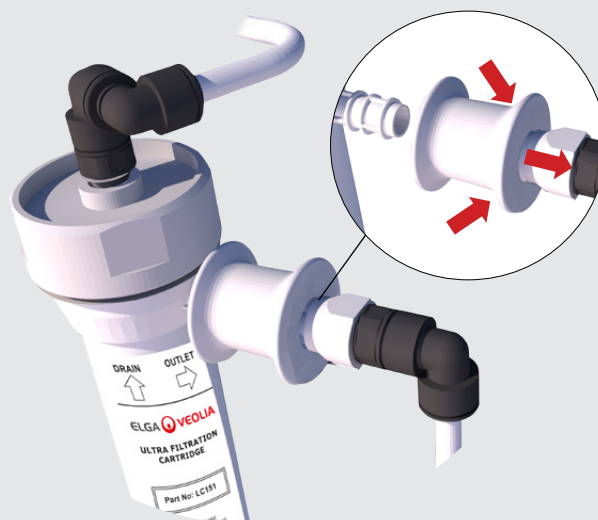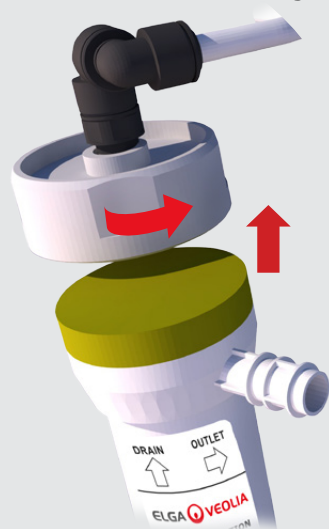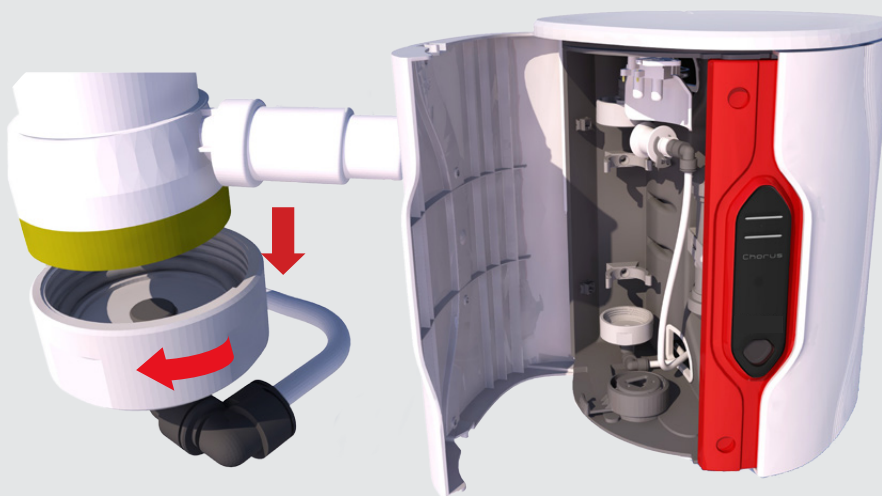**1.** Open left door, Process button off.**2.** Remove purification pack.**3.** Remove the existing LC151 from holding clips.**4.** Release the side connection, push back the sleeve of the quick connect coupling.**5.** Remove top cap, drain any liquid.**6.** Remove the bottom cap.**WARNING!**

MODULES WILL REQUIRE RINSING PRIOR TO USE.


7. Unpack new LC151 and remove transit caps .

8. Attach bottom side plug.

9. Attach the top and bottom caps (**Hand tight**).

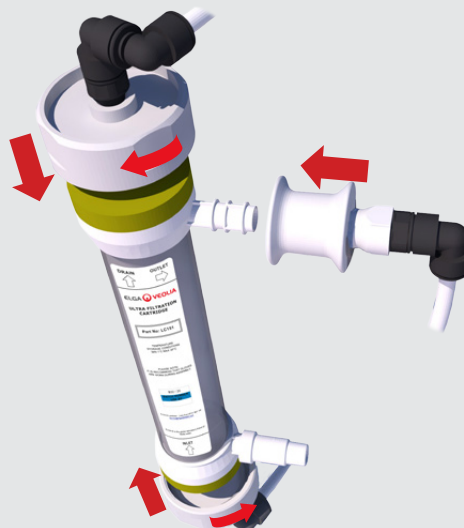10. Reconnect the quick connect coupling.

11. Install the purification pack.

12. Rinse filter. Press and hold the Process button for 15 seconds. Follow on screen instructions.




WARNING!

ULTRA FILTRATION MODULE WILL REQUIRE RINSING PRIOR TO USE.

3.5 Replacing the Ultra Mirco Filtration Module LC109

Follow the instructions below.


PC1ANRXM2

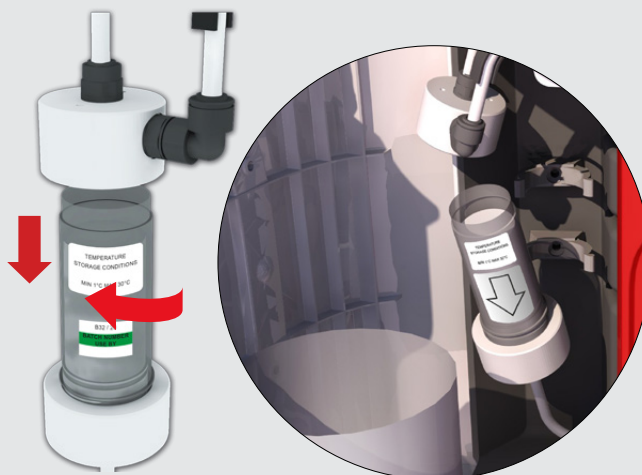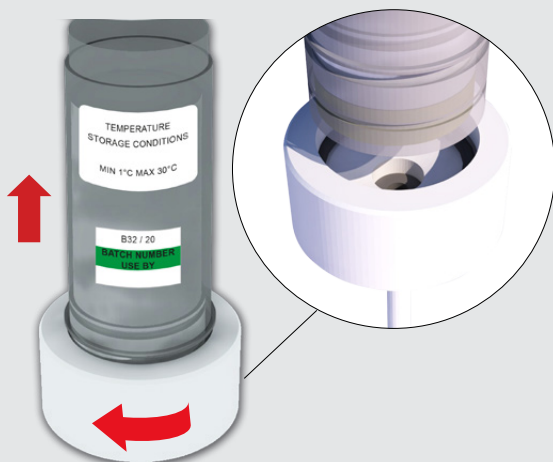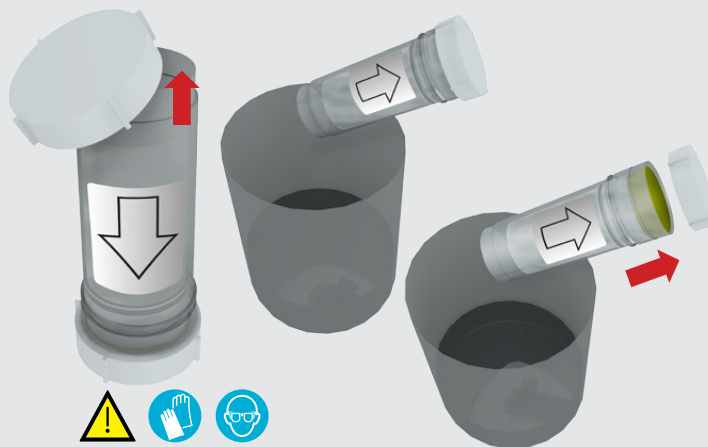**1.** Process button off, open left door.**2.** Remove purification pack.**3.** Unclip Ultra mirco filtration module LC109.**4.** Unscrew the top cap and pour the water away.**5.** Remove the bottom cap and discard LC109 filter.**6.** Unpack new LC109 Remove BOTH caps.
⚠ Caution: Pour contents to drain.**WARNING!**

THE LC109 IS SUPPLIED CONTAINING AN AQUEOUS SOLUTION OF 0.3% HYDROGEN PEROXIDE AS A BACTERIOSTAT. WEAR PROTECTIVE GLASSES AND GLOVES .


7. Attach the bottom and top caps.

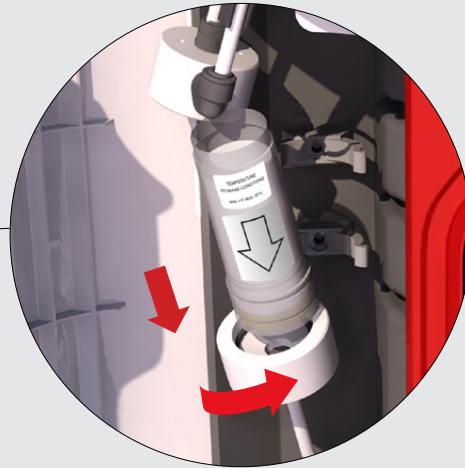Bottom cap

8. Be careful not to over tighten hand tight only.

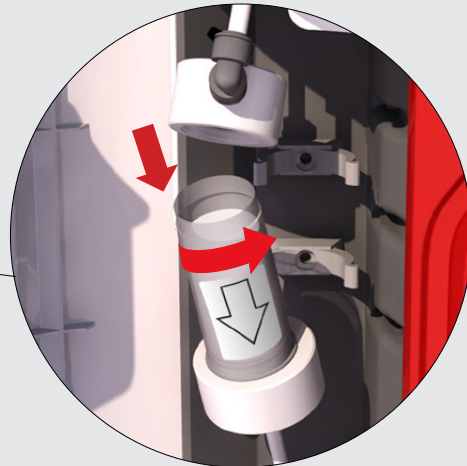Top cap

9. Insert the purification pack.

10. Rinse filter. Press and hold the Process button for 15 seconds. Follow on screen instructions.

WARNING!

ULTRA MIRCO FILTRATION MODULE WILL REQUIRE RINSING PRIOR TO USE.

3.6 Sanitisation Procedure

PURELAB® Chorus LC271 - Liquid Sanitisation - Instructions provide.

PURELAB® Chorus LC272 - Tablet Sanitisation - Instructions provide.

Please take care to ensure your order is the correct consumables items.

4 SELF HELP GUIDELINES

Display Lights

Constant	White	Signifying water purity, unit is operating correctly
Blinking	White	Consumables are nearing the end of useful life
Constant	Red	System needs immediate attention
Blinking	Red	System in clean cycle (Sanitization / Commissioning)
Blinking	Red / White	Software update

Process Button Operation

Press (At any time)	Start / Stop Modules process	Start and stops modules process
Press and hold for 15 seconds (At any time)	Enter / Exit Commissioning Procedure	To enter or exit commissioning mode / sanitization procedure. Press process button for fifteen seconds this will enter commissioning mode. Press once at anytime to pause / resume commissioning mode. This can be done using the Halo dispenser in 'Main Menu'
Press Once (At any time)	Load firmware from USB Flash devices	Automatically loads system firmware from USB flash device. Re-start on completion. This can be done using the Halo dispenser.

Default Alarm Settings

Set Temperature Alarm	Set default values from firmware file. Adjustable set point using application service module. Service module is an engineers tool, only an engineer can adjust the settings. Select to either 30°C, 35°C, 40°C. Default 35°C.
Set Water Purity Alarm Chorus 1	Set default values from firmware file. Adjustable set point using application service module. Resistivity range select from 1.0MΩ.cm, 10.0MΩ.cm or 15.0MΩ.cm. Default Alarm point 15.0MΩ.cm.
Over fill Protection Alarm	Feature activates drain valve if water level is greater than maximum and stops water entering the system to prevent a flood. When alarm is active: V2 energised. If alarm for 5mins, stop module process. Enter refill cycle on overfill alarm deactivation. Alarm status removed (Reservoir 95%)

5 HEALTH & SAFETY NOTES

WARNING!

WARNINGS ARE GIVEN WHERE FAILURE TO OBSERVE THE INSTRUCTIONS COULD RESULT IN INJURY OR FATALITY.

CAUTION!

Cautions are given where failure to observe the instructions could result in damage to the equipment, associated equipment and processes.

5.1 Environment

CAUTION!

Clean dry indoor. Temp 5-40°C. Humidity max 80% non-condensing.
Failure to follow the environmental specification could result in damage to the system.
The Chorus is not designed for use in fume cupboards where chemicals could damage the system.

WARNING!

ALL NEW UF CARTRIDGE ARE FILLED WITH A PRESERVATIVE SOLUTION TO PREVENT BACTERIAL CONTAMINATION DURING STORAGE.
THE UF CARTRIDGE WILL THEREFORE REQUIRE RINSING PRIOR TO USE.

5.2 Electricity

The appliance coupler (mains lead) or power supply connected to the rear of the unit can be removed to isolate the power supply. If access to this is restricted then it is recommended that access to the supply socket is easily available to disconnect the electrical supply.

WARNING!

ONLY USE THE APPLIANCE COUPLER (MAINS LEAD) AND POWER SUPPLY PROVIDED.
THE USE OF THESE WILL ENSURE ADEQUATE EARTH PROTECTION IS PROVIDED.

IF THE EQUIPMENT IS USED IN A MANNER NOT SPECIFIED BY ELGA, THE PROTECTION PROVIDED BY THE EQUIPMENT MAY BE IMPAIRED.

POSITION THE POWER SUPPLY SO THAT IT CANNOT COME INTO CONTACT WITH WATER

WARNING!

ALWAYS ENSURE THAT THE ELECTRICAL POWER SUPPLY IS ISOLATED BEFORE WORKING INSIDE CHORUS.

5.3 Residual Preservative

WARNING!

DURING THE COMMISSIONING CYCLE TRACE RESIDUAL PRESERVATIVE IS FLUSHED FROM THE SYSTEM.
A MATERIAL DATA SHEET IS AVAILABLE UPON REQUEST.

5.4 Ultraviolet Lamp

CAUTION!

Hold on to the UV lamp carefully to ensure it doesn't break.

CAUTION!

Take care not to touch the surface of the glass. Ideally wear gloves, handle with soft cloth and wipe the surface with alcohol before fitting into the housing.

5.5 Control of Substance Hazardous to Health (COSHH)

Material safety data sheets are available upon request.

6 TECHNICAL SPECIFICATIONS

Purified Water Specification				
Model	PURELAB® Chorus PC1LSCXM2	PURELAB® Chorus PC1ANRXM2	PURELAB® Chorus 1 PC1GSCXM2	
Delivery flow rate - maximum	2 l/m	2 l/m	2 l/m	
CO2 reduction (feed dependant)	~ 100%	~ 100%	~ 100%	
Resistivity (ionised species)	18.2 MΩ.cm	18.2 MΩ.cm	18.2 MΩ.cm	
Organics (TOC) - Typical	1 - 3 ppb	1 - 3 ppb	3 - 10 ppb	
Organics (MW >200 Dalton)	N/A	N/A	N/A	
Bacteria - Typical POU filter	<0.1 CFU/ml	<0.1 CFU/ml	<1 CFU/ml	
Bacterial Endotoxin	<0.001 EU/ml	N/A	N/A	
Particles	<0.01 um	<0.05 um	<0.2 um	
Feed Water Specification				
Source Quality	RO, DI or SDI	RO, DI or SDI	RO, DI or SDI	
Fouling Index-Maximum	1	1	1	
TDS-Maximum	N/A	N/A	N/A	
Conductivity (RO/SDI)	<30 / <1 uS/cm	<30 / <1 uS/cm	<30 / <1 uS/cm	
TOC/Organics (typical)	<50ppb	<50ppb	<50ppb	
Free/total Chlorine-Maximum	0.05ppm	0.05ppm	0.05ppm	
Chloramine - Maximum	0.05ppm	0.05ppm	0.05ppm	
Heavy Metals-Maximum (Fe/Mn)	N/A	N/A	N/A	
Silica-Maximum	2 ppm	2 ppm	2 ppm	
Temperature	1-35°C	1-35°C	1-35°C	
Flowrate (max requirement @ 15°C)	130l/h	130l/h	130l/h	
Drain requirements	2 l/m	2 l/m	2 l/m	
Maximum feed pressure	0.7 bar (10psi)	0.7 bar (10psi)	0.7 bar (10psi)	
Minimum feed pressure	0.07 bar (1psi)	0.07bar (1 psi)	0.07bar (1 psi)	

Connections

Inlet	8mm (5/16) OD tube
Outlet	8mm (5/16) OD tube
Drain	8mm (5/16) OD tube

Electrical Requirements

Model	PURELAB® Chorus PC1LSCXM2	PURELAB® Chorus PC1ANRXM2	PURELAB® Chorus 1 PC1GSCXM2
Mains Input	100-240 Vac 50-60 Hz	100-240 Vac 50-60 Hz	100-240 Vac 50-60 Hz
System control voltage	24V dc	24V dc	24V dc
Power consumption (peak demand)	60VA	60VA	60VA
Reservoir Level Connection	RJ45	RJ45	RJ45
Noise level	<45dBA	<45dBA	<45dBA

Dimensions & Weight

Model	PURELAB® Chorus PC1LSCXM2	PURELAB® Chorus PC1ANRXM2	PURELAB® Chorus 1 PC1GSCXM2
Height	435mm	435mm	435mm
Width	375mm	375mm	375mm
Depth	340mm	340mm	340mm
Supply Weight	20kg(44lb)	17kg(37lb)	17kg(37lb)
Operating Weight	22kg(49lb)	18kg(40lb)	22kg(49lb)
Installation	Bench / Wall / Floor	Bench / Wall / Floor	Bench / Wall / Floor

Dependent on feed water as part of our policy of continual improvement we reserve the right to alter the specifications given in this document.

Consumables


Part No.	Description	Typical Service Life *	Max. Shelf Life
LC210	185/254mm UV Lamp	12 - 18 Months	2 years
LC232 / LC244 / LC245 / LC246	Purification Pack	12 Months	2 years
LC109	Ultra Micro Filtration	12 Months	2 years
LC151	Ultra Filtration	12 Months	2 years
CT1	Tablets	1 every 6 months	2 years
Effersan™	Tablets	>8 Months	>8 Months

*Service life is an estimate only, and will depend on the application and feed water quality.
Please take care to ensure your order is the correct consumables items.

7 CUSTOMIZED OPERATION

It's possible to customize your Chorus by downloading additional features and changing your current settings. To find out more, please visit www.elgalabwater.com/customize

7.1 USB Data Logging To download technical data please follow instructions.

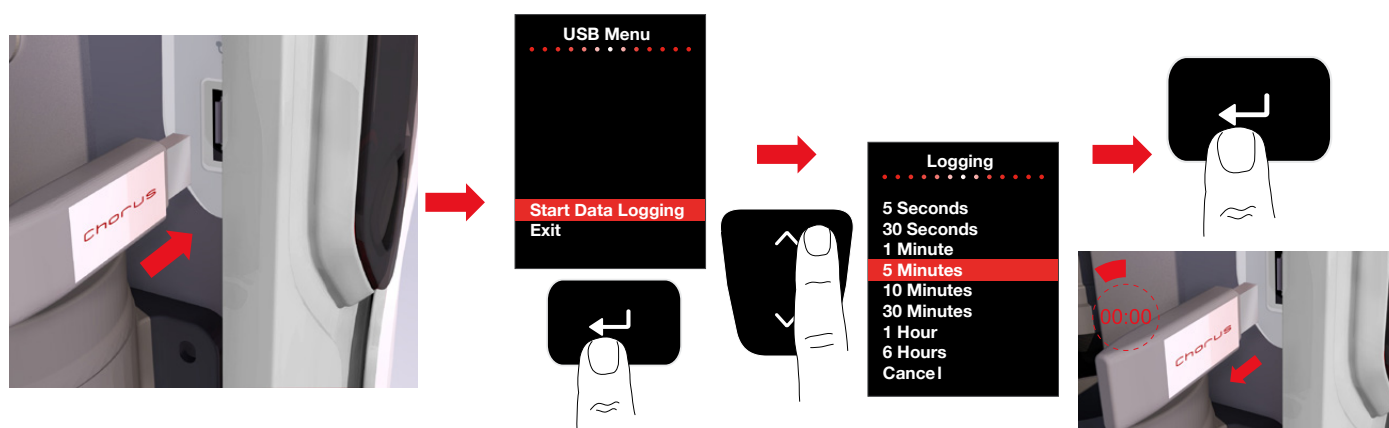ELGA® LabWater is a trading name of VWS (UK) Ltd. General Limited Warranty

VWS (UK) Ltd. warrants the products manufactured by it against defects in materials and workmanship when used in accordance with applicable instructions for a period of one year from the date of shipment for the products. VWS (UK) Ltd. MAKES NO OTHER WARRANTY, EXPRESSED OR IMPLIED. THERE IS NO WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE.

The warranty provided herein and the data, specifications and descriptions of the VWS (UK) Ltd. products appearing in VWS (UK) Ltd. published catalogues and product literature may not be altered except by express written agreement signed by an officer of VWS (UK) Ltd. Representations, oral or written, which are inconsistent with this warranty or such publications are not authorized and, if given, should not be relied upon.

In the event of a breach of the foregoing warranty, VWS (UK) Ltd. sole obligation shall be to repair or replace, at its option, any product or part thereof that proves to be defective in materials or workmanship within the warranty period, provided the customer notifies VWS (UK) Ltd. promptly of any such defect. The exclusive remedy provided herein shall not be deemed to have failed of its essential purpose so long as VWS (UK) Ltd. is willing and able to repair or replace any nonconforming VWS (UK) Ltd. product or part. VWS (UK) shall not be liable for consequential, incidental, special or any other indirect damages resulting from economic loss or property damage sustained by any customer from the use of its products.

VWS (UK) Ltd. Warranty

VWS (UK) Ltd. warrants the water systems manufactured by it, BUT EXCLUDING MEMBRANES AND PURIFICATION PACKS, against defects in materials and workmanship when used in accordance with the applicable instructions and within the operating conditions specified for the systems for a period of one year from the earlier of:

the date of installation, or the 120th day following the date of shipment.

VWS (UK) LTD. MAKES NO OTHER WARRANTY, EXPRESSED OR IMPLIED. THERE IS NO WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE. The warranty provided herein and the data, specifications and descriptions of the VWS (UK) Ltd. systems appearing in VWS (UK) Ltd. published catalogues and product literature may not be altered except by express written agreement signed by an officer VWS (UK) Ltd. Representations, oral or written, which are inconsistent with this warranty or such publications are not authorised and, if given, should not be relied upon.

In the event of a breach of the foregoing warranty, VWS (UK) Ltd. sole obligation shall be to repair or replace, at its option, any product or part thereof that proves to be defective in materials or workmanship within the warranty period, provided the customer notifies VWS (UK) Ltd. promptly of any such defect. The cost of labour for the first ninety (90) days of the above warranty period is included in the warranty; thereafter, labour cost shall be at the customer's expense. The exclusive remedy provided herein shall not be deemed to have failed of its essential purpose so long as VWS (UK) Ltd. is willing and able to repair or replace any nonconforming VWS (UK) Ltd. system or component part. VWS (UK) Ltd. shall not be liable for consequential, incidental, special or any other indirect damages resulting from economic loss or property damage sustained by any customer from the use of its process systems.

Products or components manufactured by companies other than VWS (UK) Ltd. or its affiliates

("Non - VWS (UK) Ltd. products") are covered by the warranty, if any, extended by the Product manufacturer. VWS (UK) Ltd. hereby assigns to the purchaser any such warranty; however VWS (UK) LTD. EXPRESSLY DISCLAIMS ANY WARRANTY WHETHER EXPRESSED OR IMPLIED, THAT THE NON - VWS (UK) Ltd. PRODUCTS ARE MERCHANTABILITY OR FIT FOR A PARTICULAR PURPOSE.

NOTICE

VWS (UK) Ltd. is constantly striving to improve its products and services. Consequently, the information in this document is subject to change without notice and should not be construed as a commitment by VWS (UK) Ltd.

Also VWS (UK) Ltd. assumes no responsibility for any errors that may appear in this document. This manual is believed to be complete and accurate at the time of publication. In no event shall VWS (UK) Ltd. be liable for incidental or consequential damages in connection with or arising from the use of this manual.

VWS (UK) Ltd. warrants its products against defects in materials and workmanship as described in the Warranty statement on the preceding pages.

The Labwater Specialists

ELGA is the global laboratory water brand name of Veolia. The information contained in this document is the property of VWS (UK) Ltd, trading as ELGA LabWater, and is supplied without liability for errors or omissions. © VWS (UK) Ltd.
2017 - All right reserved. ELGA®, PURELAB®, ADEPT® Labpure® and PureSure® are registered trademarks of VWS (UK) Ltd.

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© 2017 ELGA Labwater / VWS (UK) Ltd. All rights reserved.